## Bio-Link – BLX 使用手冊

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- *Crosslinking of DNA and RNA by covalently binding nucleic acids to nitrocellulose or nylon membranes for Southern, Northern, dot or slot blots procedures (1-2-3)*

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- *Nicking of Ethidium Bromide stained DNA in agarose gel (4)*
- *Partial restriction endonuclease digestion by formation of thymine dimers (5)*
- *RecA mutation screening (6)*
- *Elimination of PCR contamination (7)*
- *The BIO-LINK® BLX is useful for other applications such as UV sterilization, UV curing of polymers, etc...*

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## →保固 Warranty

Our products (except Compact Flash®, light tubes and filters) are warranted against faulty construction or defective material for a period of TWO YEARS from the date of supply. Our products are not warranted for damage due to carelessness, incorrect use or bad maintenance.

The following defects are also specifically excluded:

- defects caused by improper operation

Repair or modification done by anyone other than Vilber Lourmat or an authorized agent.

-corrosion caused by improper solvents or samples.

-use of spare parts supplied by anyone other than Vilber Lourmat.

-damage caused by accident or misuse.

-damage caused by disaster.

This instrument should not be modified or altered in any way. Modification or alteration of this instrument will:

1. void the manufacturer's warranty.
2. void the conformity certifications.
3. create a potential safety hazard.

The Compact Flash®, the tubes and the filters are not covered by our warranty. The use of consumable products or non-original spare parts not recommended by our service department is at the user's own risk and therefore automatically invalidates the warranty. Tubes, filters, batteries and consumable products are not included in the warranty. We reserve the right to decide where the faulty goods will be repaired (in our workshop or elsewhere), and whether or not the faulty part is to be replaced; all other freight charges

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incurred being at the cost of the purchaser. Returned goods will not be accepted for repair unless previous written authorisation is obtained from our service department. A request for authorisation must be accompanied by an itemised list of products, model numbers and the corresponding invoice numbers under which they were originally shipped. All returned goods should have a certificate of decontamination. The Buyer must bear all costs and risks incurred during the transportation of the goods from their collection at Vilber Lourmat warehouse. In the case VILBER LOURMAT incorporates some devices or equipment from another supplier in the manufacture of its products, the extent and the duration of the warranty will be those conceded by the suppliers or sellers. Manufacturer cannot be held responsible for any loss, bodily injury or material accident incurred by any failure of this supply, whatever the origin of this failure may be. The responsibility of Manufacturer is strictly limited to its staff and to its own supplies. In the case of dispute, only the commercial court of Meaux (France) shall be competent, even in third party claims proceedings or when there are several co-defendants. NOTE: Vilber Lourmat is not responsible for any injury or damage caused by use of this instrument for purposes other than those for which it is intended, or by modifications of the instrument not performed by Vilber Lourmat.

#### Decontamination, collection and elimination of waste for France only

The buyer ensures and finances the decontamination, the collection and the disposal of waste electrical and electronic equipment (WEEE) under the conditions provided in the Articles 21 and 22 of the Decree No. 2005-829 dated of 20 July 2005. In France, for tubes recycling, contact the company Recylum, [www.recylum.com](http://www.recylum.com) Improper disposal may be harmful to the environment and human health.

## → Declaration of conformity

	<p>The materials comply with the requirements of the EC Directive 89/336/EEC, 73/23/EEC and EN 61010-1 (electro-magnetic compatibility and low voltage).</p> <p>The electro-magnetic susceptibility has been chosen at a level that gains proper operation in residential areas, on business and light industrial premises and on small-scale enterprises, inside as well as outside of the buildings. All places of operation are characterized by their connection to the public low voltage power supply system.</p>
---	---

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